

ZB-5MSplus Drugs of AbuseColumn: Zebtron[™] ZB-5MSplus, GC Cap. Column 30 m x 0.25 mm x 0.25 μ m, Ea

Phase: Proprietary Pesticides Phase

Dimensions: 30 meters x 0,25 mm x 0,25 μ m

Order No: 7HG-G030-11

Oven Profile: 70 for 1 min to 150 at 30 °C/min to 320 at 10 for 5.5 min

Carrier Gas: Constant Flow Helium, 1,4 mL/min


Injection: Splitless :1 1 μ L @ 250°C

Detection: Mass Spectrometer (MS) (250°C)

Detector Info: 70 for 1 min to 150 at 30 °C/min to 320 at 10 for 5.5 min

Analyst Note: MSD: 45 -450 da
Sample @ 25ppm

Products used in this application:

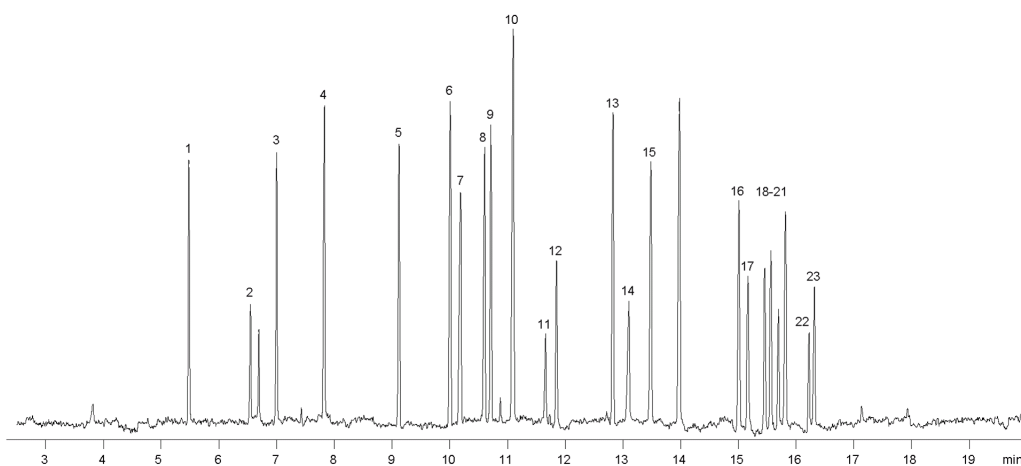

ZB-5MSplus Drugs of Abuse

ANALYTES:

- 1 Nicotine
- 2 MDA
- 3 Mdma
- 4 MBDB
- 5 Meperidine
- 6 Benzamphetamine
- 7 Glutethimide
- 8 PCP
- 9 Doxylamine
- 10 Tramadol
- 11 Procaine
- 12 Methadone
- 13 Topiramate
- 14 Cocaine
- 15 Cocaethylene
- 16 CBD
- 17 Codeine
- 18 Ethylmorphine
- 19 Diazepam
- 20 Hydrocodone
- 21 D9-THC
- 22 6-mam
- 23 Oxycodone

